Package 'licoread'

May 8, 2025

Title Reads Raw Files from Li-COR Gas Analyzers Version 0.2.0 Date 2025-05-08 Description Reads raw files from Li-COR gas analyzers and produces a dataframe that can directly be used with 'fluxible' <https://www.com/actionality.com/actional //cran.r-project.org/package=fluxible>. **License** GPL (≥ 3) **Encoding** UTF-8 RoxygenNote 7.3.2 **Suggests** knitr, rmarkdown, testthat (>= 3.0.0) Config/testthat/edition 3 Imports dplyr, jsonlite, lubridate, purrr, readr, rlang, stringr, tibble, tidyr, yaml URL https://jogaudard.github.io/licoread/ BugReports https://github.com/jogaudard/licoread/issues VignetteBuilder knitr NeedsCompilation no Author Joseph Gaudard [aut, cre] (ORCID: <https://orcid.org/0000-0002-6989-7624>) Maintainer Joseph Gaudard <joseph.gaudard@pm.me> **Depends** R (>= 4.1.0) **Repository** CRAN

Date/Publication 2025-05-08 16:50:09 UTC

Contents

data_82z																													2
data_name_	82z																												3
fluxible_81x	κ	•	•	•	•	•	•	 •	•	•	•	•	•		•	•	•	•	•	•	•				•	•	•	•	3

fluxible_82z	4
licoread	4
licoread_auto	5
licoread_to_fluxible	6
list_gases	7
metadata_82z	. 7
names_df	. 8
oneobs_81x	8
oneobs_82z	9
read_81x_onefile	. 9
units_82z	10
	11

Index

data_82z

to read the raw data

Description

to read the raw data

Usage

```
data_82z(filepath, data_file, data_name, gases, filename)
```

Arguments

filepath	name and path to the 82z archive
data_file	name of the file with raw data
data_name	vector of colnames
gases	list of gases
filename	name of the 82z archive

Value

a long df with the actual data contained in the data file

data_name_82z create colnames for data tibble

Description

create colnames for data tibble

Usage

```
data_name_82z(filepath, data_file)
```

Arguments

filepath	name and path to the 82z archive
data_file	name of the file with raw data

Value

a character vector or the colnames of the data file

fluxible_81x makes df from 81x files compatible with fluxible

Description

makes df from 81x files compatible with fluxible

Usage

fluxible_81x(df, focus_gas, id_cols, datetime_col)

Arguments

df	input dataframe from licoread
focus_gas	gas to select
id_cols	columns to identify unique fluxes
datetime_col	column containing datetime information

Value

a df with the focus gas column renamed as "f_conc" and f_fluxid in chronological order of datetime

fluxible_82z

Description

makes df from 82z files compatible with fluxible

Usage

```
fluxible_82z(df, focus_gas)
```

Arguments

df	input dataframe from licoread
focus_gas	gas to select

Value

an unnested df with only the selected gas

licoread

reads Li-COR files in a given location

Description

reads Li-COR files in a given location

Usage

```
licoread(
  location,
  file_type = "auto",
  file_type_list = c("82z", "81x", "auto"),
  data_file = "data.csv",
  meta_file = "metadata.json",
  regex_file = "(\\w*-)*\\w*(?=([.]82z$))",
  sample = FALSE
)
```

licoread_auto

Arguments

location	location of the files
file_type	type of file (82z or 81x). If "auto" (default), the function will try to detect it by itself.
file_type_list	list of file types
data_file	name of the file with raw data
<pre>meta_file</pre>	name of the file with meta data
regex_file	regex expression matching the name of the 82z file. Here in case the user has a different than the default and for easier updates.
sample	sample = n randomly selects n files to be imported. This allows for testing the setup before importing a potentially large list of files which will take time and be difficult to handle.

Value

a tibble (nested or not depending on raw data) containing all the data from the raw files present at the location provided

Examples

```
path_82z <- system.file("extdata/82z", package = "licoread")
licoread(path_82z)</pre>
```

licoread_auto finds out the file type for licoread

Description

finds out the file type for licoread

Usage

```
licoread_auto(file_list)
```

Arguments

file_list list of files found in the location

Value

a single character string indicating the file type

licoread_to_fluxible selects the focus gas and makes a df to use in fluxible

Description

selects the focus gas and checks the columns needed for the fluxible workflow

Usage

```
licoread_to_fluxible(
  df,
  focus_gas,
  datetime_col,
  id_cols = c("File Name", "Obs#"),
  file_type = "auto",
  file_type_list = c("82z", "81x", "auto")
)
```

Arguments

df	input dataframe from licoread
focus_gas	gas to select
datetime_col	column containing datetime information if date and time are in two different columns, provide a character vector of the form c("date", "time")
id_cols	columns to identify unique fluxes
file_type	type of file (82z or 81x). If "auto" (default), the function will try to detect it by itself.
file_type_list	list of file types

Value

an unnested df with only the selected gas

Examples

```
path_82z <- system.file("extdata/82z", package = "licoread")
gas_df_82z <- licoread(path_82z)
licoread_to_fluxible(gas_df_82z, "LI-7810_CH4_DRY",
datetime_col = c("LI-8250_DATE", "LI-8250_TIME"))</pre>
```

list_gases

Description

prints a character vector of the gases present in the dataset. This is to be used to obtain the exact names of the gases before using licoread_to_fluxible.

Usage

```
list_gases(df, file_type = "auto", file_type_list = c("82z", "81x", "auto"))
```

Arguments

df	the dataframe as imported with licoread
file_type	type of file (currently works only for 82z). If "auto" (default), the function will try to detect it by itself.
file_type_list	list of file types

Value

a character vector of the gases present in the dataset

Examples

```
path_82z <- system.file("extdata/82z", package = "licoread")
gas_df_82z <- licoread(path_82z)
list_gases(gas_df_82z)</pre>
```

metadata_82z read meta data file inside 82z archive

Description

read meta data file inside 82z archive

Usage

```
metadata_82z(filepath, meta_file)
```

Arguments

filepath	name and path to the 82z archive
<pre>meta_file</pre>	name of the file with meta data

Value

a tibble with the metadata from one observation

names_df

Description

to get a vector of names of a df, matching a regex

Usage

```
names_df(df, regname = "name\\d")
```

Arguments

df	the df to get the names from
regname	the regex expression to match

Value

a df with the names of the meta df

oneobs_81x reading a single measurement from 81x file

Description

reads a single measurement from a licor .81x file

Usage

```
oneobs_81x(start, end, all_obs, file)
```

Arguments

start	line number at which the measurement starts
end	line number at which the measurement ends
all_obs	list of all the lines from the full file
file	filepath to the 81x files

Value

a df with 1 row with the meta data of the measurement and raw data nested

oneobs_82z

Description

to read one measurement from the 82z archive

Usage

oneobs_82z(filepath, data_file, meta_file, regex_file)

Arguments

filepath	path to the 82z archive
data_file	name of the file with raw data
<pre>meta_file</pre>	name of the file with meta data
<pre>regex_file</pre>	regex expression matching the name of the 82z file. Here in case the user has a different than the default and for easier updates.

Value

a tibble with all the data and metadata from one observation (one file)

read_81x_onefile reads 81x licor file

Description

reads a .81x file with several measurements

Usage

```
read_81x_onefile(file)
```

Arguments

file filepath the the 81x file to read

Value

a nested tibble with the meta data from each measurements as row and the raw data nested

units_82z

Description

to create a nested tibble with the units of data

Usage

units_82z(filepath, data_file, data_name, filename)

Arguments

filepath	name and path to the 82z archive
data_file	name of the file with raw data
data_name	vector of colnames
filename	name of the 82z archive

Value

a tibble with the units of the variables contained in the raw data

Index

data_82z, 2
data_name_82z, 3
fluxible, 6
fluxible_81x, 3
fluxible_82z, 4
licoread, 3, 4, 4, 6, 7
licoread_auto, 5
licoread_to_fluxible, 6, 7
list_gases, 7
metadata_82z, 7
names_df, 8

oneobs_81x, 8 oneobs_82z, 9

read_81x_onefile,9

units_82z, 10